

SULFIDE TEST PROCEDURE



The Sulfide test procedure utilizes eXact® Strip Micro and eXact® reagent to accurately achieve a Sulfide reading with the eXact iDip® Smart Photometer System®. For tips on achieving best accuracy, and additional information regarding features of the eXact iDip® photometer and the eXact iDip® app, please refer to the eXact iDip® Smart Guide. You can download a copy on our website at exactidip.com. **Test procedures can vary from test to test. Read full instructions and watch instructional video within the app.**


YOU WILL NEED

- eXact iDip® photometer
- Your smartphone/tablet with the eXact iDip® app installed and Sulfide S2 purchased in the Store
- 1 eXact® Strip Micro S (Part No. 486818-A)
- 4 drops eXact® Reagent S (Part No. 486818-B)
- Water sample for rinsing and testing

1 SELECT CUSTOMER

- Select **'Customers'** from the Home screen.
- Choose customer from list
 - OR** Tap **'Add customer from contacts'** and choose customer from your contacts list
 - OR** Tap **'+'** to create a new customer
- Android users: If no address is found, tap "No addresses found"
- Verify customer has been selected

Note: To take full advantage of the GPS and Data Storage features, each test result must be linked to a contact. Individual users may select their contact listing.

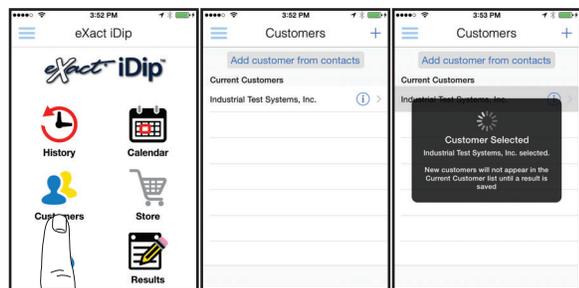

a

b


c

2 POWER ON EXACT IDIP®

Press the **ZERO/ON** button to power on the eXact iDip® photometer.

3 SELECT BLUETOOTH TEST


Tap the menu slide out **'≡'** and select **'Bluetooth Test'** from the choices shown.

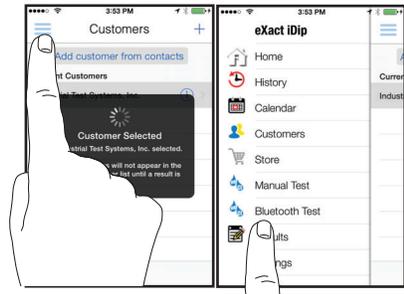

4 CONNECT EXACT IDIP®


The eXact iDip® app will automatically connect to the most recently used eXact iDip® photometer. If not, select your eXact iDip® from the bottom of the screen.

Note: Always connect your eXact iDip® photometer via the Bluetooth® connection within the app. If you experience an issue connecting your device, check that your smartphone/tablet's Bluetooth® is turned on and your device is compatible.

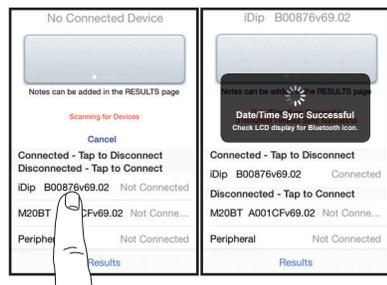

5 SELECT TEST

Tap **'SELECT'** Test' at the top and select **Sulfide S2** (if this is not visible, purchase the test from the Store). The eXact iDip® photometer and eXact iDip® app will both display the test selection for Sulfide.

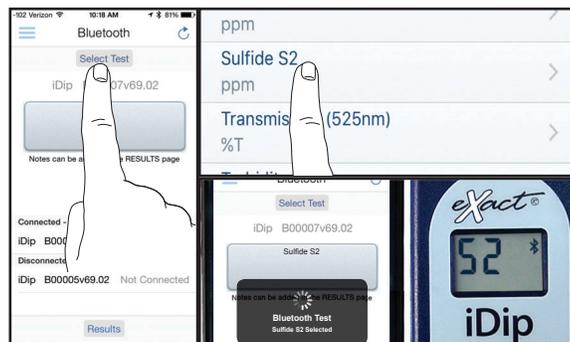

6 FILL CELL

Before testing, rinse CELL and clean with brush thoroughly. Finally, rinse the cell 3 times with the water sample to be tested, then **FILL** cell to capacity to begin test (refer to video in app).



7 ADD DROPS

Add **4 drops** of eXact® Reagent S (Part No. 486818-B). Ensure the reagent bottle remains vertical while adding drops.




8 COVER & MIX

Place the Cell Cover onto the CELL. Turn the meter **upside-down** and shake the meter up and down three times to mix the sample.



9 ZERO METER

Press **ZERO/ON** and the eXact iDip® photometer display reads 0 PPM indicating the meter is ready for testing.

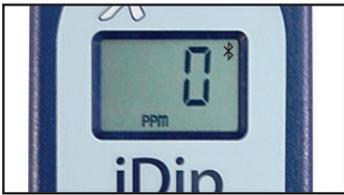

10 REMOVE STRIP

Remove one eXact® Strip Micro S (Part No. 486818) and set in a dry, convenient place. Immediately replace cap on bottle.



11 DIP STRIP


Press **READ** to initiate a 20 second countdown and simultaneously **DIP** the eXact® strip in the sample, gently touching the bottom of the cell. Use a gentle constant back and forth motion (2 strokes per second) until the timer displays "1". Remove and discard the strip.

The eXact iDip® photometer will now display a 180 second count-up timer. Make sure your smart device does not go to sleep.



12 CAP CELL AND READ RESULTS

Place the Cell Cover onto the CELL and **READ** result displayed on the eXact iDip® photometer and in the eXact iDip® app.

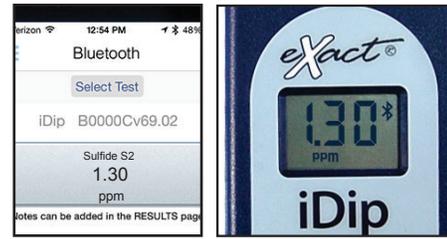

13 CLEAN CELL WITH BRUSH

After testing, rinse CELL immediately and clean with brush to remove reagents which coat the CELL wall.



TIPS

- Due to the manufacturing process, you may find one or two strips that are noticeably smaller or larger in width than the normal strips in the bottle. These should be discarded. Using these strips may give unreliable results.
- Before running test unlock Sulfide test by purchasing in the Store.
- After the 20 second countdown, the display will immediately start counting up from 1 to 40.
- The calibration of the meter is based on the water sample temperature above 20°C (68°F). If the water sample is below 20°C (68°F), the strip has to dip in the sample for an additional 10 seconds.
- To obtain optimal accuracy when testing in direct sunlight, use the Cell Cover when zeroing and reading the sample.
- For results as Hydrogen Sulfide (H₂S), multiply the resulting value by 1.06.
- It is recommended to use the Cell Cleaning Brush with Distilled White Vinegar, 0.1N HCl, or Muriatic Acid to clean the CELL after each test.
- Do not run this test before or after Total Iron test.
- Be careful while dipping the strip (step 11). Do not spill the sample from the CELL.

MANAGE YOUR RESULTS

Save, send, and share results instantly using the eXact iDip® app. For step-by-step instructions, refer to the eXact iDip® Smart Guide. You can download a pdf copy on-line at exactidip.com.



eXact iDip

